




# Handabfüllstation für flüssigen Stickstoff

Zur Benutzung dieser Abfüllstation müssen die folgenden Voraussetzungen erfüllt sein:

1. Gültiges FS-Sicherheitstraining (DOOR) „Tiefkalte Gase“
2. Persönliche Schutzausrüstung (Kryohandschuhe und Gesichtsschutzschild)
3. Einweisung durch den jeweiligen Hallenmeister
4. Anwesenheitspflicht während des Vorgangs
5. Die Betriebsanweisung für tiefkalte Gase ist zu beachten
6. Kleine Dewars in den Edelstahltopf stellen

Es kann einige Minuten dauern, bis Flüssigkeit aus dem goldenen Korken austritt.

Es darf keinerlei Flüssigkeit auf den Boden tropfen. Die Epoxydbeschichtung platzt sofort.

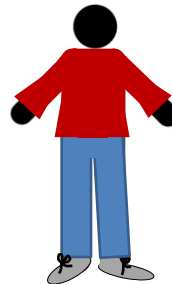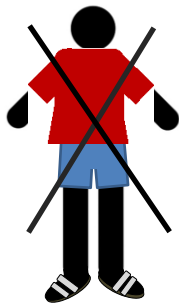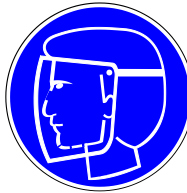

# Filling Station for Liquid Nitrogen

In order to use this filling station, the following preconditions must be fulfilled:

1. Valid safety training (DOOR) „Cryogenic Gases“
2. Personal protective equipment (cryo gloves and face shield)
3. Instruction by the responsible hall supervisor
4. Compulsory attendance during filling
5. The operating instructions cryogenic liquids must be followed
6. Small dewars must be placed inside the large vessel

It might take several minutes before liquid nitrogen is flowing out of the „golden cork“. Be patient.

It has to be avoided that liquid droplets touch the floor.

The epoxy coating will be destroyed immediately.

